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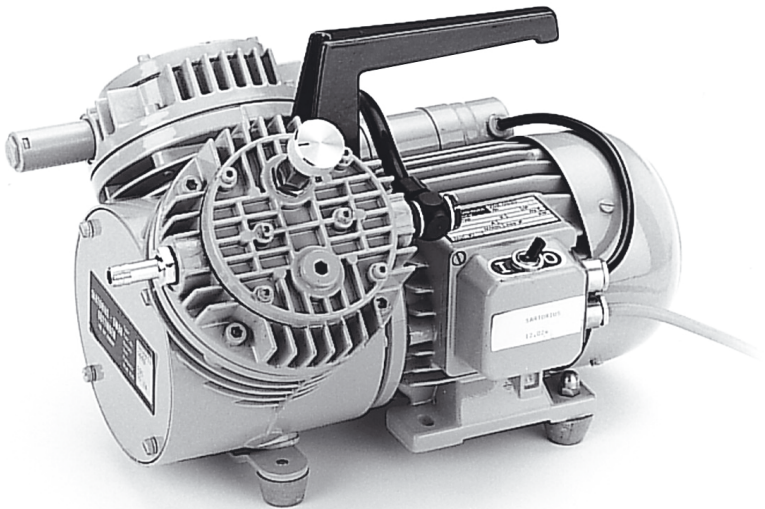
## Contents

<b>1 Specifications</b> .....	4
<b>2 Important Notes</b> .....	4
<b>3 Maintenance</b> .....	5
3.1 Changing the Diaphragm.....	5
3.2 Changing the Reed Valves.....	6
<b>4 Recommended Accessories</b> .....	7
<b>5 Spare Parts</b> .....	7
<b>EC/EU Declaration of Conformity</b> .....	8
<b>FCC Supplier's Declaration of Conformity</b> .....	9

The Sartorius Stedim Biotech Electric Diaphragm Vacuum Pump is especially convenient for using multibranch vacuum manifolds. Built to last, it runs quietly while providing a high performance. The pump comes equipped with a carrying handle; rubber feet; built-in power supply with line cord (2 m); muffler; serially connected twin pump heads on the suction side; and hose connector.

The reducing valve allows the vacuum to be regulated. The final vacuum is 13 mbar (10 torr). The pump's maximum suction rate is 26 l/min.

The internationally recognized protection rating IP 44\* for the pump motor conforms to the German standard VDE 0530, Section 5. The electrical design of the pump motor complies with the German standard VDE 0530, Section 1.

\* described in the IEC Publication 144 and entails protection against contact and splashing

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16615: 110 V, 60 Hz

Diaphragm: neoprene

Max. pumping rate (at atmospheric pressure): 26 l/min.

Final vacuum: 13 mbar (10 torr)

Motor output: 120 W

Power consumption: 1.8 amps

Max. permissible ambient temperature: 40°C

Weight: approx 9.8 kg

Dimensions (mm): 338 x 250 x 225

Connections: connector for 9 mm vacuum hose

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## 2 Important Notes

1. Before you plug in the pump, check the manufacturer's label to make sure the voltage rating matches your local line voltage (mains).
2. The ambient temperature may not exceed 40°C. Pumps installed in equipment by the customer must have sufficient ventilation inside and outside the cabinet.
3. Do not start the pump under load (vacuum).
4. The pump may only be used for air, gases and non-aggressive vapors, but not liquids.
5. Protect the pump against jets and gushes of water and against excessive dust.

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## 3 Maintenance

The Sartorius Stedim Biotech Diaphragm Vacuum Pump does not require regular maintenance. The ball bearings are oversized and sufficiently lubricated with a special grease to last for the pump's entire life. The diaphragms and valves are the only parts subject to wear. However, changing these parts does not require any special technical expertise and can be easily done on-site within a few minutes.

### 3.1 Changing the Diaphragm (see Fig.1).

3.1.1 Remove the connecting tube between the diaphragm heads (1) by loosening the threaded nut (4) and removing the tube (3) from one side. Alternatively, you can remove the tube completely with the screws from the diaphragm head (unscrew).

3.1.2 Use a pencil to mark the position of the diaphragm head (1) in relation to the housing (11).

3.1.3 Untighten the 4 Allen screws (12), and remove the diaphragm head.

3.1.4 Now loosen the countersunk screw, and remove the retainer plate, star lock washer and diaphragm one by one.

3.1.5 Remove the end cover plate (6) for the motor fan from the right side of the pump (remove the 4 cheese head screws) and turn the disk flywheel until the connecting rod is in mid-position. Next, install the new diaphragm.

3.1.6 Now replace the retainer plate first with the flat side facedown on the diaphragm followed by the star lock washer, making sure that it is properly aligned. Carefully but securely tighten the countersunk screw.

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3.1.7 Next, replace the diaphragm head in the marked position and uniformly tighten the Allen screws in opposite diagonal pairs.

3.1.8 Check that the pump runs smoothly by rotating the disk flywheel one complete turn and then replace the end cover plate.

3.1.9 Reattach the tube to connect both diaphragm heads.

## **3.2 Changing the Reed Valves**

### **I. Lower Valve**

3.2.1 Remove the diaphragm head as described in steps 3.1.1 to 3.1.3.

3.2.2 Loosen the cheese head screw, and exchange the spring steel reed valves.

### **II. Upper Reed Valve**

3.2.3 Loosen the 6 Allen screws (10) and remove the cover plate (9) and gasket. If necessary, remove the paint from the edge of the cover beforehand (use a screw-driver to scrape it off).

3.2.4 Loosen the cheese head screw, and exchange the reed valves.

3.2.5 Retighten the cheese head screw, and replace the cover plate with a new gasket. Retighten the 6 Allen screws. Remount the diaphragm head, and check that the pump runs smoothly as described in 3.1.7 and 3.1.8.

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16610 Woulff'sche Flasche, 500 ml  
16623 Gummivakuumschlauch (1 m)

## 4 Recommended Accessories

16610 Woulff's bottle, 500 ml  
16623 Rubber vacuum hose (1 m)

## 5 Ersatzteile

6986017

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- 2 Membranen
- 4 Ventildfedern
- 2 Kopfdichtungen

## 5 Spare Parts

6986017

Spare parts kit consisting of:

- 2 diaphragms
- 4 reed valves
- 2 gaskets for the diaphragm heads

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*The person authorised to compile the technical file:*

Sartorius Lab Instruments GmbH & Co. KG  
International Certification Management  
37070 Goettingen, Germany

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Sartorius Lab Instruments GmbH & Co. KG  
Goettingen, 2016-12-20

Dr. Reinhard Baumfalk  
Vice President R&D

Dr. Dieter Klausgrete  
Head of International Certification Management

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*This declaration certifies conformity with the above mentioned EC and EU Directives, but does not guarantee product attributes. Unauthorised product modifications make this declaration invalid. The safety information in the associated product documentation must be observed.*

# FCC Supplier's Declaration of Conformity

## FCC Supplier's Declaration of Conformity

**Device type** Pump

**Model** 16615, 16695, 16694-1-60-06, 16694-1-60-22

**Party issuing Supplier's Declaration of Conformity /  
Responsible Party – U.S. Contact Information**

Sartorius Corporation  
5 Orville Dr Suite 200  
11716 Bohemia, NY  
USA  
Telephone: +1.631.254.4249

**FCC Compliance Statement**

This device complies with Part 15 of the FCC Rules. Operation is subject to the following two conditions: (1) This device may not cause harmful interference, and (2) this device must accept any interference received, including interference that may cause undesired operation.

**Information to the user**

Note: This equipment has been tested and found to comply with the limits for a **class B** digital device, pursuant to part 15 of the FCC Rules. These limits are designed to provide reasonable protection against harmful interference in a residential installation. This equipment generates, uses and can radiate radio frequency energy and if not installed and used in accordance with the instructions, may cause harmful interference to radio communications. However, there is no guarantee that interference will not occur in a particular installation. If this equipment does cause harmful interference to radio or television reception, which can be determined by turning the equipment off and on, the user is encouraged to try to correct the interference by one or more of the following measures:

- Reorient or relocate the receiving antenna.
- Increase the separation between the equipment and receiver.
- Connect the equipment into an outlet on a circuit different from that to which the receiver is connected.
- Consult the dealer or an experienced radio/TV technician for help.

Connections between the device and peripherals must be made using shielded cables in order to maintain compliance with FCC radio frequency emission limits.

Any modifications made to this device that are not approved by Sartorius may void the authority granted to the user by the FCC to operate this equipment.

Sartorius Lab Instruments GmbH & Co. KG  
Otto-Brenner-Strasse 20  
37079 Goettingen, Germany

Phone: +49.551.308.0  
www.sartorius.com

The information and figures contained in these instructions correspond to the version date specified below.

Sartorius reserves the right to make changes to the technology, features, specifications and design of the equipment without notice. Masculine or feminine forms are used to facilitate legibility in these instructions and always simultaneously denote the other gender as well.

Copyright notice:

This instruction manual, including all of its components, is protected by copyright. Any use beyond the limits of the copyright law is not permitted without our approval. This applies in particular to reprinting, translation and editing irrespective of the type of media used.

© Sartorius Germany

Last updated:  
10 | 2018