



Virus Counter[®] 3100 Instrument

The Solution for Rapid, Direct, Biologically Relevant Virus Quantification

Features

- **Fast results**
Near real-time quantification of viruses for highly efficient manufacturing optimization
- **Biologically relevant readout**
Specific, direct, high affinity labeling systems
- **Versatility**
Broadly reactive reagents for use with a number of viruses
- **Ease of use**
Software-assisted system operation and data analysis (21 CFR Part 11 capabilities)



Product Information

The Virus Counter[®] 3100 platform is a powerful analytical tool that provides near real-time virus counts. The system can be used throughout virus growth and purification processes in a number of applications. The platform enables users to identify optimal harvesting times, validate processing steps, and define final formulations with speed and confidence.

Viruses are first labeled with specific reagents, either using a direct dye approach or high affinity monoclonal antibodies. The samples are then subjected to precise and intense laser illumination. This approach allows rapid, direct and specific quantification of total virus particles in a sample.

The different detection methods of the Virus Counter[®] platform offer the right solution for individual quantification needs.

Description

The Virus Counter[®] 3100 platform, that includes the instrument, software, consumables and reagents, was purpose-built to enumerate viruses and other small particles. In this system, viruses are first directly labeled with fluorescent dyes or fluorescently-labeled antibody-based reagents. Within the instrument, virus samples are guided into a fluidic stream and focused to a small core of moving fluid. A laser is used to excite the fluorophores associated with the viruses as they pass through the laser area. The Virus Counter[®] instrument detects the emitted light and uses the signals together with the measured sample flow rate to calculate total particle concentration.

Applications

- Optimization of viral growth conditions to increase product yield
- Real-time insights into virus titers in bioreactors for the determination of optimal harvesting times
- Optimization of processing steps and tracking of virus titers through downstream purification steps
- More accurate characterization of final vector product for increased safety and efficiency
- Critical benefits in a number of industries:
 - Vaccine production
 - Protein expression
 - Viral therapeutics in gene and cell therapy
 - Antiviral development

System Operation

The operation of the Virus Counter[®] 3100 instrument is software guided. For instrument performance verification and sample analysis several consumables are used. Two are directed specifically to verify instrument performance while three are washing solutions that ensure minimal crossover between samples. It is critical to use the individual consumables for instrument maintenance and washing steps to achieve optimal performance. All consumables are color coded for easy discrimination

Fluidics

Accurate quantification requires that the virus sample stream passes through a laser interrogation point and that the particles traverse that point temporally separated. Differential flow rates of the sheath and sample lead to accelerated laminar flow as the core stream passes in front of the laser spot and collection lens. This hydrodynamic focusing of the sample stream ensures that sample particles travel through the interrogation point at the peak of laser intensity.

Optics

As the laser illuminates the narrow sample stream within the flow cell, microscope optics are used to collect fluorescence from stained particles as they pass through the interrogation point. A dichroic mirror is used to separate the two fluorescence signals. Bandpass filters spectrally filter and isolate each channel's fluorescence. The fluorescence emissions are detected by photomultiplier tubes (PMTs).

Virus Signal Detection

Signals collected from the PMTs are processed in real time to convert the observed fluorescence information into a quantitative measure of virus particles per ml (vp/ml). For Combo Dye[®] reagent, a mixture of a protein and a nucleic acid dye, light of two different wavelengths is emitted upon laser excitation. Both detection channels record fluorescence events that exceed a predetermined amplitude threshold that identifies events above background. While all events are collected, only the events that occur simultaneously in both channels are used to indicate that a virus particle was present in the light path. Therefore, only simultaneous events for Combo Dye[®] reagent are included in the concentration determination. ViroTag[®] results represent fluorescence emission from a fluorochrome-conjugated antibody reagent in one channel. Instrument quantification limits are 5×10^5 to 1×10^9 virus particles/ml.

Data Analysis

Experimental data is automatically collected and stored within a SQL database installed on the Virus Counter[®] laptop. This data can be viewed and analyzed using the Virus Counter[®] Software data analysis feature. Alternatively, data can be exported as a .csv file and viewed using third party data analysis packages.

Service

Reliability and smooth operation of equipment are ensured through proper installation | configuration and regular maintenance by Sartorius Service experts.



Technical Data

Laser Product Classification	Class 1: Complies with EN 60825-1 and US 21 CFR 1040.10 Note: Internally, an enclosed Class 3B laser diode module is used per EN 60825-1.
Instrument Air Source	Industrial Grade Compressed Gas – Clean Dry Air (CDA) or Nitrogen: 12-15 psi (103 kPa)
Operating Environment	+18°C – +27°C (+65°F – +80°F), < 60% Relative Humidity
Indoor Outdoor Use Category	Indoor Use Only
Weight	13.5 kg, 29.8 lbs. (dry weight)
Dimensions (W × D × H)	43 cm × 43 cm × 28 cm (17" × 17" × 11")
Power Requirements – Computer	19.5 Vdc 2.3 A
Power Requirements – Instrument	12 Vdc 5 A
External Power Supply – Computer	Input: 100 – 240 V (+/-10%) ~50 – 60 Hz 150 VA, Inlet Plug IEC320/C6, mating cord IEC320/C5
External Power Supply – Instrument	Input: 100 – 240 V (+/-10%) ~50 – 60 Hz 1.5 VA, Inlet Plug IEC320/C14, mating cord IEC320/C13 Output: 12 Vdc 5 A
Data I/O Communications Port	USB 2.0
Ambient Pollution Degree	Pollution Degree 2 per EN 61010-1
Noise Level	<70 dB (A)
Enclosure Protection Class	Ordinary Protection Only (Not protected against harmful Ingress of Moisture). IP20, NEMA-1

Ordering Information

Instrument	Order Number
Virus Counter [®] 3100 Instrument	VIR-92341

Virus Counter Reagents

Reagent	Target	Order Number	Quantity
ViroTag [®] AAV2-3	Intact AAV2 and AAV3 capsids	VIR-92117	200 samples per kit
ViroTag [®] BCBV	Budded Baculovirus	VIR-92108	200 samples per kit
ViroTag [®] INVA	Influenza A Viruses (H1 and H3 subtype)	VIR-91151	200 samples per kit
ViroTag [®] INVB	Influenza B viruses	VIR-91152	200 samples per kit
ViroTag [®] VSVG	VSV-G Protein	VIR-92332	200 samples per kit
Combo Dye [®]	Universal stain for enveloped viruses	VIR-92333	200 samples per kit

Virus Counter Consumables

Consumable	Order Number	Quantity
VC3100 Sheath Fluid	VIR-92302	1 L bottle
VC Wash Fluid	VIR-92296	1 L bottle
VC Sample Dilution Buffer	VIR-92298	1 bottle
VC Combo Dye [®] Dilution Buffer	VIR-92305	1 bottle

Sales and Service Contacts

For further contacts, visit www.sartorius-stedim.com

Europe

Germany

Sartorius Stedim Biotech GmbH
August-Spindler-Strasse 11
37079 Goettingen

Phone +49.551.308.0

Sartorius Stedim Systems GmbH
Robert-Bosch-Strasse 5-7
34302 Guxhagen

Phone +49.5665.407.0

France

Sartorius Stedim FMT S.A.S.
ZI des Paluds
Avenue de Jouques - CS 91051
13781 Aubagne Cedex

Phone +33.442.845600

Sartorius Stedim France SAS
ZI des Paluds
Avenue de Jouques - CS 71058
13781 Aubagne Cedex

Phone +33.442.845600

Austria

Sartorius Stedim Austria GmbH
Modectcenterstrasse 22
1030 Vienna

Phone +43.1.7965763.18

Belgium

Sartorius Stedim Belgium N.V.
Rue Colonel Bourg 105
1030 Bruxelles

Phone +32.2.756.06.80

Hungary

Sartorius Stedim Hungária Kft.
Kagyló u. 5
2092 Budakeszi

Phone +36.23.457.227

Italy

Sartorius Stedim Italy S.r.l.
Via dell'Antella, 76/A
50012 Antella-Bagno a Ripoli (FI)

Phone +39.055.63.40.41

Netherlands

Sartorius Stedim Netherlands B.V.

Phone +31.30.60.25.080

filtratie.nederland@sartorius-stedim.com

Poland

Sartorius Stedim Poland Sp. z o.o.
ul. Wrzesinska 70
62-025 Kostrzyn

Phone +48.61.647.38.40

Russian Federation

LLC "Sartorius Stedim RUS"
Vasilyevsky Island
5th line 70, Lit. A
199178 St. Petersburg

Phone +7.812.327.53.27

Spain

Sartorius Stedim Spain, S.A.U.
Avda. de la Industria, 32
Edificio PAYMA
28108 Alcobendas (Madrid)

Phone +34.913.586.098

Switzerland

Sartorius Stedim Switzerland AG
Ringstrasse 24 a
8317 Tagelswangen

Phone +41.52.354.36.36

U.K.

Sartorius Stedim UK Ltd.
Longmead Business Centre
Blenheim Road, Epsom
Surrey KT19 9 QQ

Phone +44.1372.737159

Ukraine

LLC "Sartorius Stedim RUS"
Post Box 440 "B"
01001 Kiev, Ukraine

Phone +380.44.411.4918

Africa

Republic of South Africa

Sartorius South Africa (Pty) Ltd
Unit 4, Alphen Square South
853 16th Road
Midrand 1685, RSA

Phone +27.11.315.5444

Americas

USA

Sartorius Stedim North America Inc.
565 Johnson Avenue
Bohemia, NY 11716

Toll-Free +1.800.368.7178

Argentina

Sartorius Argentina S.A.
Int. A. Ávalos 4251
B1605ECS Munro
Buenos Aires

Phone +54.11.4721.0505

Brazil

Sartorius do Brasil Ltda
Avenida Senador Vergueiro 2962
São Bernardo do Campo
CEP 09600-000 - SP- Brasil

Phone +55.11.4362.8900

Mexico

Sartorius de México, S.A. de C.V.
Libramiento Norte de Tepotzotlan s/n,
Colonia Barrio Tlacateco,
Municipio de Tepotzotlan,
Estado de México,
C.P. 54605

Phone +52.55.5562.1102

leadsmex@sartorius.com

Asia | Pacific

Australia

Sartorius Stedim Australia Pty. Ltd.
Unit 5, 7-11 Rodeo Drive
Dandenong South Vic 3175

Phone +61.3.8762.1800

China

Sartorius Stedim (Shanghai)
Trading Co., Ltd.
3rd Floor, North Wing, Tower 1
No. 4560 Jinke Road
Zhangjiang Hi-Tech Park
Pudong District
Shanghai 201210, P.R. China

Phone +86.21.6878.2300

Sartorius Stedim (Shanghai)
Trading Co., Ltd.
Beijing Branch Office
No. 33 Yu'an Road
Airport Industrial Park Zone B
Shunyi District, Beijing 101300

Phone +86.10.8042.6501

Sartorius Stedim (Shanghai)
Trading Co., Ltd.
Guangzhou Branch Office
Room 1105
Xing Guang Ying Jing Building
No. 119, Shui Yin Road
Yue Xiu District, Guangzhou 510075

Phone +86.20.3836.4193

India

Sartorius Stedim India Pvt. Ltd.
#69/2-69/3, NH 48, Jakkasandra
Nelamangala Tq
562 123 Bangalore, India

Phone +91.80.4350.5250

Japan

Sartorius Stedim Japan K.K.
4th Fl., Daiwa Shinagawa North Bldg.
8-11, Kita-Shinagawa 1-chome
Shinagawa-ku, Tokyo, 140-0001 Japan

Phone +81.3.4331.4300

Malaysia

Sartorius Stedim Malaysia Sdn. Bhd.
Lot L3-E-3B, Enterprise 4
Technology Park Malaysia
Bukit Jalil
57000 Kuala Lumpur, Malaysia

Phone +60.3.8996.0622

Singapore

Sartorius Stedim Singapore Pte. Ltd.
10 Science Park Rd
The Alpha #02-13/14
Singapore Science Park II
Singapore 117684

Phone +65.6872.3966

South Korea

Sartorius Korea Biotech Co., Ltd.
8th Floor, Solid Space B/D,
PanGyoYeok-Ro 220, BunDang-Gu
SeongNam-Si, GyeongGi-Do, 463-400

Phone +82.31.622.5700



◀ www.sartorius-stedim.com