

Instructions for Use

Microsart[®] ATMP Extraction

Bacterial, fungi, and yeasts DNA extraction kit

Prod. No. SMB95-2001

Reagents for 50 extractions

For use in research and quality control

Manufactured by:



Minerva Biolabs GmbH
Schkopauer Ring 13
12681 Berlin
Germany

Symbols

LOT

Lot No.

REF

Order No.



Expiry date



Store at



Contains reagents for
50 extractions



Manufacturer

Contents

1. Intended Use	5
2. Explanation of the Test	5
3. Test Principle	5
4. Notes on the Test Procedure	6
4.1 Handling and Equipment Recommendations . . .	7
5. Reagents	8
6. Needed but not Included	9
7. Specimen	12
Sample Collection and Storage	12
8. Test Procedure	13
8.1 Recommendation for product release testing . . .	13
8.2 DNA Extraction Process	13
9. Short Instructions	14
10. Related Products	16



1. Intended Use

Microsart® ATMP Extraction kit introduces a state-of-the-art DNA extraction method for DNA extraction from cell cultures and cell culture derived biologicals, like autologous transplants and other advanced therapy medicinal products (ATMP), for subsequent DNA amplification via PCR.

2. Explanation of the Test

To achieve highest sensitivity and to avoid inhibitory effects in PCR testing, a DNA extraction is recommended. For most test materials, DNA extraction methods are available providing templates suitable for PCR. However, most of the DNA extraction kits available on the market are not free of DNA contaminations. Microsart® ATMP Extraction introduces a unique DNA extraction method, which eliminates the risk of DNA contaminations, facilitating the detection of bacteria, fungi, and yeasts in cell culture and ATMPs via PCR.

The extraction procedure can be performed within 1 hour. In contrast to the culture method, samples do not need to contain vital material, as all intact particles (e.g. live, dormant, non-culturable etc.) are detected.

3. Test Principle

Microsart® ATMP Extraction kit was optimized for the extraction of genomic bacterial or fungal DNA from different sample matrices including cell culture samples. The contamination risk has been reduced to a minimum due to less handling steps.

An internal amplification control DNA from the Microsart® ATMP Bacteria or Microsart® ATMP Fungi kit can be added to the sample prior DNA extraction to monitor the extraction process by detecting false negative results which can occur due to improper DNA extraction or PCR inhibition. Alternatively, the internal control DNA can be added directly to the Mastermix during PCR setup.

4. Notes on the Test Procedure

1. For in vitro use in research and quality control. This kit may be disposed of according to local regulations.
2. This kit should be used only by trained persons. You should wear a clean lab coat and use disposable gloves at all times while performing the assay.
3. To avoid DNA cross-contaminations, the complete test must be performed under sterile and DNA-free conditions (see chapter 4.1 for detailed information).
4. In case of working with living strains, the local regulatory requirements for S2 labs must be considered.
5. Attention: by aliquoting and repeatedly freezing and thawing your samples, you run a high risk of contamination. This should therefore be avoided if possible.
6. This extraction kit has been validated with 1 ml starting volume. If you use less than 1 ml it must be ensured that 99 cfu can be detected in the appropriate volume.
7. This kit is not validated for the extraction of mycoplasma DNA.
8. This leaflet must be widely understood for a successful use of Microsart® ATMP Extraction kit. The reagents supplied should not be mixed with reagents from different lots but used as an integral unit. The reagents of the kit should not be used beyond their shelf life.
9. Any deviation from the test method can affect the results.
10. For each test setup, at least one negative extraction control should be included. Positive controls facilitate the evaluation of the test.
11. The controls should be carried out in the same manner as the samples.
12. Participation in external quality control programs, such as those offered by Minerva Biolabs GmbH (www.minerva-biolabs.com), is recommended.

4.1 Handling and Equipment Recommendations

To avoid false positive results due to improper handling the following actions are recommended:

1. To perform the test under sterile and DNA-free conditions, we recommend the use of an isolator/glovebox with an airlock.
2. The isolator/glovebox should be cleaned thoroughly with PCR Clean™ (Minerva Biolabs, Prod. No. 15-2025) or PCR Clean™ Wipes (Minerva Biolabs, Prod. No. 15-2001) before and during the working process.
3. All materials which are introduced into the isolator/glovebox should be cleaned thoroughly with PCR Clean™. Don't forget to clean the airlock with PCR Clean™. Pipettes and gloves should be cleaned thoroughly with PCR Clean™ Wipes prior and during the process.
4. Avoid working above open tubes and avoid air turbulences due to rapid movements.
5. Be careful when opening the tubes. Do not touch the inner surface of the lid.

5. Reagents

Each kit contains reagents for 50 extractions. The expiry date of the unopened package is marked on the package label. The kit components are stored at ambient temperature until use.

Kit Component Label Information	50 Extractions Order No. SMB95-2001	Cap Color
Lysis Buffer	2 x 13 ml	transparent
Suspension Buffer	4 x 1.5 ml	violet
Processing Tubes	50 Tubes	

The lot specific Certificate of Analysis can be downloaded from the manufacturer's website (www.minerva-biolabs.com).

6. Needed but not Included

Microsart® ATMP Extraction kit contains reagents for sample collection and DNA extraction. General industrial supplies and reagents, usually available in PCR laboratories are not included:

Consumables

- Laboratory gloves
- PCR Clean™ (Minerva Biolabs, Prod. No. 15-2025) and PCR Clean™ wipes (Minerva Biolabs, Prod. No. 15-2001)
- DNA-free pipette filter tips (we recommend Biosphere® filter tips from Sarstedt: 0.5-20 µl, Prod. No. 70.1116.210; 2-100 µl, Prod. No. 70.760.212; 20-300 µl, Prod. No. 70.765.210; 100-1000 µl, Prod. No. 70.762.211)
- DNA-free PCR reaction tubes (PCR 8-SoftStrips with attached caps from Biozym are recommended: 0.1 ml Low Profile, Prod. No. 710975 and 0.2 ml High Profile, Prod. No. 710970)

Equipment

- Isolator/glovebox (further information, supplier and prices are available on request, please contact PCR@sartorius.com)
- Heat block with optional shaking function
- Microcentrifuge for 1.5 ml reaction tubes (Centrisart A-14, Prod. No. A-14-1EU)
- Vortex Mixer
- Pipettes
 - mechanical
 - 0.5 – 10 µl Sartorius Prod. No. LH-729020
 - 10 – 100 µl Sartorius Prod. No. LH-729050
 - 100 – 1000 µl Sartorius Prod. No. LH-729070
 - or electrical
 - 0.2 – 10 µl Sartorius Prod. No. 735021
 - 10 – 300 µl Sartorius Prod. No. 735061
 - 50 – 1000 µl Sartorius Prod. No. 735081
- Rack for 1.5 ml tubes


For subsequent PCR analysis, the following consumables and equipment are required additionally:

- DNA PCR detection system for the detection of bacteria: we recommend the Microsart® ATMP Bacteria kit (Sartorius Prod. No. SMB95-1008), or the Microsart® RESEARCH Bacteria kit (Sartorius Prod. No. SMB95-1009/1010).
- DNA PCR detection system for the detection of fungi and yeasts we recommend the Microsart® ATMP Fungi kit (Sartorius Prod. No. SMB95-1012), or the Microsart® RESEARCH Fungi kit (Sartorius Prod. No. SMB95-1013/1014).

All these kits are qPCR-based methods, designed for the detection from cell cultures and cell culture derived biologicals.

- qPCR device with filter sets for the detection of the fluorescence dyes FAM™ and ROX™ and suitable for 25 µl PCR reaction volumes
- DNA-free PCR reaction tubes that must be free from DNA for the specific qPCR device (PCR 8-SoftStrips with attached caps from Biozym are recommended: 0.1 ml Low Profile, Prod. No. 710975 and 0.2 ml High Profile, Prod. No. 710970)
- Minicentrifuge for PCR-tubes
- Isolator/glovebox (for PCR-setup)
- Vortex mixer
- Set of 3 pipettes
- Rack for 1.5 ml tubes
- Rack for PCR tubes

Schematical overview of technical setup and experimental design:

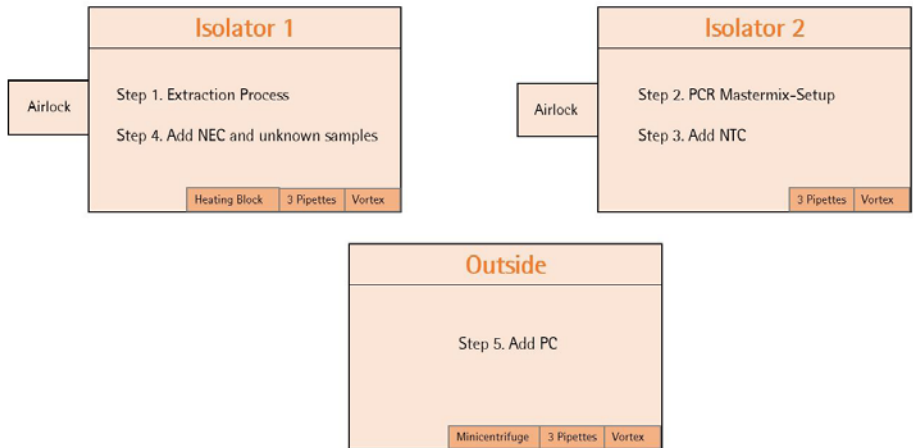


It is also possible to connect Isolator 1 and Isolator 2 via an airlock so that you can transfer the PCR tubes after Step 3 directly from Isolator 2 into Isolator 1. Please note that in this case you would need an additional airlock for Isolator 2.

7. Specimen


Sample Collection and Storage

The kit has been validated using a maximum cell concentration of 10^6 cells/ml. Notably, the assay can be performed with different types of cell culture-derived material. Therefore, the optimal sampling parameters, like volume or cell number, can vary according to the specific characteristics of the sample (e.g. medium, cell type) and may require optimization of the procedure.

1. max. 1 ml of cell culture or cell culture supernatant liquid material is transferred into a provided DNA-free 1.5 ml processing tube (transparent cap).
2. Spin down for 15 minutes at a speed of at least 16,200 x g to sediment particles. Attention: Make sure to position the tubes in the centrifuge in order to obtain a pellet on the back side of the tube, as described in the figure below.
3. Discard the supernatant carefully and completely as described in the figure below. Proceed to DNA extraction. If DNA extraction cannot be performed immediately, freeze samples at ≤ -18 °C. Repeated freezing and thawing should be avoided. Attention: Samples can only be inactivated or frozen after this sample collection step.

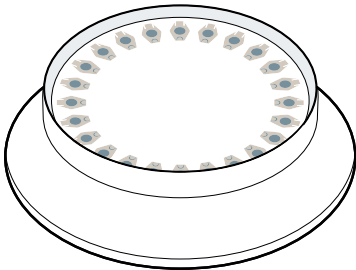

Make sure to position the tubes with the back side toward the outside of the rotor in order to obtain a pellet on the back wall of the tube.

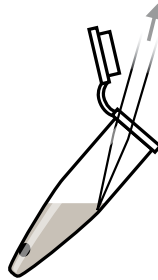

Slowly discard all the supernatant without disturbing the pellet.

8. Test Procedure

8.1 Recommendation for product release testing

The extraction process should be carried out with a negative extraction control (NEC) and samples in duplicates (= 3 extractions for 1 product).

8.2 DNA Extraction Process

-
1. Add 500 µl Lysis Buffer (transparent cap) to cell pellet.
Recommended for users of Microsart® ATMP Bacteria and Microsart® ATMP Fungi detection kit: The Internal Control DNA which is included in the detection kit can also be used to monitor the extraction process. Add 20 µl Internal Control DNA to the sample, vortex briefly and proceed with step 2 as described. No additional Internal Control DNA is required for the PCR reaction mix.
 2. Vortex vigorously for at least 30 seconds until pellet is completely dissolved.
 3. Heat at 80°C (optional with shaking) for 10 minutes.
 4. Spin down at 16,200 x g for 10 minutes.
Attention: make sure to position the tubes in the rotor as indicated on the figure in chapter 7.
 5. Remove supernatant carefully and completely following the explanations in paragraph 7. Make sure not to disturb or aspirate the pellet in the process.
Attention: There is a higher risk of inhibition in PCR analysis if residues remain in the tube.
 6. Add 100 µl Suspension Buffer (violet cap) and dissolve the DNA by thorough vortexing.
-

Extracts can be stored for 6 days at +2 to +8 °C. If long term storage is required, store at ≤ -18 °C. Repeated freezing and thawing should be avoided.

9. Short Instructions

1. Sample Collection

+ 1 ml sample material
processing tube

☉ 15 min ≥ 16,200 xg

discard supernatant

store at ≤ -18 °C
or
proceed to DNA Extraction

2. DNA Extraction

+ 500 µl Lysis Buffer (transparent cap)
optional: add Internal Control DNA from Microsart® ATMP Bacteria or Microsart® ATMP Fungi

⚡ ≥ 30 sec vigorously
🕒 80 °C, 10 min
☉ 16,200 g, 10 min

remove supernatant carefully

+ 100 µl Suspension Buffer (violet cap)
⚡ ≥ 30 sec vigorously

proceed to PCR

🕒 incubate
⚡ vortex
☉ centrifuge
+ add

storage at ambient temperature
after rehydration ≤ -18 °C

Appendix

Limited Product Warranty

This warranty limits our liability for replacement of this product. No warranties of any kind, express or implied, including, without limitation, implied warranties of merchantability or fitness for a particular purpose, are provided. Minerva Biolabs shall have no liability for any direct, indirect, consequential, or incidental damages arising out of the use, the results of use, or the inability to use this product.

Trademarks

Microsart is a registered trademark of Sartorius Stedim Biotech, GmbH. PCR Clean is a trademark of Minerva Biolabs, GmbH.

Last technical revision: 2019-11-15



10. Related Products

Detection Kits for qPCR

SMB95-1001/1002	Microsart® AMP Mycoplasma	25/100 tests
SMB95-1003/1004	Microsart® ATMP Mycoplasma	25/100 tests
SMB95-1005/1006	Microsart® RESEARCH Mycoplasma	25/100 tests
SMB95-1009	Microsart® RESEARCH Bacteria	25 tests
SMB95-1008	Microsart® ATMP Bacteria	100 tests
SMB95-1007	Microsart® ATMP Sterile Release	10 samples
SMB95-1012	Microsart® ATMP Fungi	100 tests
SMB95-1014/1013	Microsart® RESEARCH Fungi	25/100 tests

Microsart® Calibration Reagent, 1 vial, 10⁸ genomes / vial (bacteria, including Mollicutes)

SMB95-2021	Mycoplasma arginini
SMB95-2022	Mycoplasma orale
SMB95-2023	Mycoplasma gallisepticum
SMB95-2024	Mycoplasma pneumoniae
SMB95-2025	Mycoplasma synoviae
SMB95-2026	Mycoplasma fermentans
SMB95-2027	Mycoplasma hyorhinis
SMB95-2028	Acholeplasma laidlawii
SMB95-2029	Spiroplasma citri
SMB95-2030	Bacillus subtilis
SMB95-2031	Pseudomonas aeruginosa
SMB95-2032	Kocuria rhizophila
SMB95-2033	Clostridium sporogenes
SMB95-2034	Bacteroides vulgatus
SMB95-2035	Staphylococcus aureus
SMB95-2036	Mycoplasma salivarium

Microsart® Calibration Reagent, 1 vial, 10⁶ genomes / vial (fungi)

SMB95-2044	Candida albicans
SMB95-2045	Aspergillus brasiliensis
SMB95-2046	Aspergillus fumigatus
SMB95-2047	Penicillium chrysogenum
SMB95-2048	Candida glabrata
SMB95-2049	Candida krusei
SMB95-2050	Candida tropicalis

Microsart® Validation Standard, 3 vials each, 10 CFU / vial (Mollicutes)

SMB95-2011	Mycoplasma arginini
SMB95-2012	Mycoplasma orale
SMB95-2013	Mycoplasma gallisepticum

SMB95-2014	Mycoplasma pneumoniae
SMB95-2015	Mycoplasma synoviae
SMB95-2016	Mycoplasma fermentans
SMB95-2017	Mycoplasma hyorhinis
SMB95-2018	Acholeplasma laidlawii
SMB95-2019	Spiroplasma citri
SMB95-2020	Mycoplasma salivarium

Microsart® Validation Standard, 6 vials each, 99 CFU / vial (bacteria* and fungi)

SMB95-2005	Bacillus subtilis
SMB95-2006	Pseudomonas aeruginosa
SMB95-2007	Kocuria rhizophila
SMB95-2008	Clostridium sporogenes
SMB95-2009	Bacteroides vulgatus
SMB95-2010	Staphylococcus aureus
SMB95-2037	Candida albicans
SMB95-2038	Aspergillus brasiliensis
SMB95-2039	Aspergillus fumigatus
SMB95-2040	Penicillium chrysogenum
SMB95-2041	Candida glabrata
SMB95-2042	Candida krusei
SMB95-2043	Candida tropicalis

* except for Mollicutes

DNA Extraction Kit

SMB95-2003	Microsart® AMP Extraction (for mycoplasma)	50 extractions
56-0002	Proteinase K**	50 extractions

PCR Clean™ **

15-2025	DNA Decontamination Reagent, spray bottle	250 ml
15-2200	DNA Decontamination Reagent, refill bottles	4x 500 ml

PCR Clean™ Wipes**

15-2001	DNA Decontamination Wipes	50 wipes
15-2002	DNA Decontamination Wipes, refill sachets	5 x 50 wipes

** Distributed by Minerva Biolabs

Sartorius Stedim Biotech GmbH
August-Spindler-Str. 11
37079 Goettingen, Germany

Phone +49.551.308.0
Fax +49.551.308.3289
www.sartorius-stedim.com

Copyright by
Sartorius Stedim Biotech GmbH,
Goettingen, Germany.
All rights reserved. No part of
this publication may be reprinted
or translated in any form or by any
means without the prior written
permission
of Sartorius Stedim Biotech GmbH.
The status of the information,
specifications and illustrations
in this manual is indicated by
the date given below.
Sartorius Stedim Biotech GmbH
reserves the right to make changes
to the technology, features,
specifications and design of the
equipment without notice.

Status:
November 2019,
Sartorius Stedim Biotech
GmbH, Goettingen, Germany

Printed in Germany on paper that
has been bleached without any use
of chlorine. | W
Material No.: 1000054395
Ver.11 | 2019